

LDMS Specimen Management Codes and Volume for MTN 003

Use the table below as a guide when entering 003 specimens in the LDMS. For each test, enter the corresponding codes and volume (primary and aliquot) in the LDMS Specimen Management module.

| Test | Primary | Primary Additive | Primary Volume | Primary Units | Aliquot Derivative | Aliquot Sub Add/Der | Aliquot Volume | Aliquot Units |
|---------------------|---------|------------------|----------------|---------------|--------------------|---------------------|----------------|---------------|
| Vaginal Swabs | VAG | PBS | 2 | EA | SWB | N/A | 1 | EA |
| Endocervical Swabs | CXS | PBS | 2 | EA | CXS | N/A | 1 | EA |
| Vaginal Gram Stains | VAG | NON | 2 | EA | GRS | GRS | 1 | EA |
| Plasma for storage | BLD | EDT | Variable | ml | PL 1/2 | N/A | 1.0–2.0 | ml |