

# NICHD LDMS Reference Guide

## Labels Module: Verifying Labels

Once labels are generated and printed, verify that they have been printed successfully.

This can be completed by opening Notepad (or similar) and scanning a barcode. If the label was printed successfully, the global specimen ID from the barcode will appear in Notepad.

It is expected to see the “FSQ” prefix included to the global specimen ID. When generating 2D support LDMS barcodes, LDMS will append the “FSQ” prefix to the LDMS global specimen ID.

FSQ9999-0002SC00-001



ABCDEFGH  
9999-0002SC00-001  
123ABC456 STUDYABC  
04/Aug/2021 08:30  
BLD EDT PL2 N/A  
1.00 ML 2 Vst